



Home for Cells and Beyond

PGmatrix-BioInk-hiPSC Using Guide

PGmatrix-BioInk for hiPSC (PGbioInk) takes the advantage of its unique features including shear-thinning, rapid gel-recovery (self-healing), and tunable gel strength capability, offering the best cyto-compatible hydrogel to meet advanced tissue engineering and organ biofabrication via 3D bioprinting techniques without using light or chemical crosslinking agent. PGbioInk Kit consists of a vial of PGmatrix-BioInk nanofiber solution, a vial of PGworks trigger solution, and a vial of PGgrow-hiPSC solution. PGbioInk kit creates a microenvironment mimicking in vivo ECM accordingly for cells growing and migrating in 3D manner. All operating procedures can be completed at room temperature or 37°C in neutral pH condition.

NOTE 1: PGmatrix-Spheroid (PG-S) kit with similar mechanical properties and cyto-compatibility as the PGbioink kit can also be used for 3D bioprinting following the PGbioInk-hiPSC protocol.

PRODUCT: PepGel™ PGmatrix-BioInk for hiPSC Research Kit
CONTENT: PGmatrix-BioInk solution, PGworks solution, and PGgrow-hiPSC
QUANTITY: 20 mL of 4% PGmatrix-BioInk, 4 mL of PGworks, and 500 µL PGgrow-hiPSC
10 mL of 4% PGmatrix-BioInk, 2 mL of PGworks, and 250 µL PGgrow-hiPSC
6 mL of 4% PGmatrix-BioInk, 1 mL of PGworks, and 150 µL PGgrow-hiPSC

STORAGE: Stored at 4°C

FOR IN VITRO RESEARCH USE ONLY. PLEASE READ MATERIAL USING AGREEMENT FOR MORE DETAILS. FOR IN VIVO TEST, PLEASE ASK FOR PG IN VIVO PRODUCTS.

CONTACT: PepGel LLC
101 BIVAP Bldg.
1980 Kimball Ave
Manhattan, KS 66506
E-mail: info@pepgel.com
Phone: 785-226-6917
Fax: 785-532-7193
Web: www.pepgel.com

To Order: customerservice@pepgel.com, or online www.pepgel.com

PGmatrix-BioInk for hiPSCs Using Protocol

1. Cell suspension preparation: To prepare cell medium stock solution, thaw PGgrow and add into E8 complete medium at ratio 1:1000 v/v (PGgrow : E8 complete medium) and use within two weeks after dilution. Suspend cells/spheroids in stock medium to prepare cell suspension. Recommend final cell seeding density in bioink hydrogel for printing is 2×10^5 - 4×10^5 cell/mL (6000-7000 spheroids/mL). For example: to prepare 2% PGbioInk hydrogel with 4×10^5 cell/mL, cell density of cell suspension needs to be 9.5×10^5 cell/mL.

2. Mixing ratios: The PGbioInk solution contains 4% W/V standard peptides, specially designed for hiPSCs. **Tables 1** presents mixing ratios of PGbioink solution and PGworks to obtain final gel concentration of 0.5%, 1%, 2%, and 3% as reference, respectively. Then follow procedure #3 below for bioink preparation.

Table 1: Examples of Mixing ratios of 4% PGbioInk solution, PGworks* and cell suspension (PGworks is always 16% of the total volume of PGbioInk solution.)**

Final concentration***	PGbioInk solution (μL)	PGworks (μL)	Cell suspension	Total volume (μL)
3%	750	120	130	1000
2%	500	80	420	1000
1%	250	40	710	1000
0.5%	125	20	855	1000

NOTES: * add PGworks to your cell suspension **FIRST** before you mix PGbioInk solution with cell suspension.

** Complete E8 medium with PGgrow is recommended for 3D culturing of hiPSC in bioprinted constructs.

***Bioink from 0.5 – 1% peptide concentration is suitable for ink-jet printing or manually pipetting for small construct desirable for Lab-on-Chip research, and bioink from 1.5% above is for extrusion printer or other printing devices.


3. Protocol for PGbioInk for hiPSC Preparation

1. Bring the PGbioInk solution and PGworks solution to room temperature (15 - 25 °C) or 37 °C (37 °C water bath)


PepGel

Home for Cells and Beyond

2. Suspend cells in completed E8 medium with PGgrow, then add PGworks solution to the cell suspension according to the Mixing Ratio in **Table 1**, pipet well without introducing air bubbles (always immersing pipet tip in cell solution during pipetting).

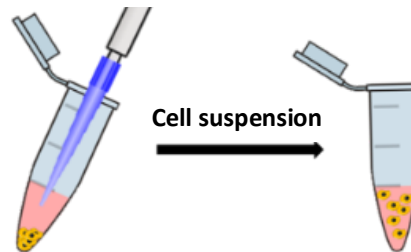

3. Mix the PGbioInk solution carefully into the cell suspension from step 2 at the Mixing Ratio following **Table 1** (pipet well without introducing air bubbles). Transfer the mixture into the center of the bioink container (i.e., syringe or pipet).
4. Incubate the bioink container at 37°C (5% CO₂) for 30 min to complete the gelation, then it is ready for bioprinting. (**NOTE:** for hydrogel infusion purpose, the mixture from step 3 can be used, and then perform step 4 for gelation after infusion).
5. Culture the bioprinted construct at 37°C (5% CO₂) using complete E8 medium supplemented with PGgrow. Replace medium to feed cells depends on cell growth without disturbing the construct.

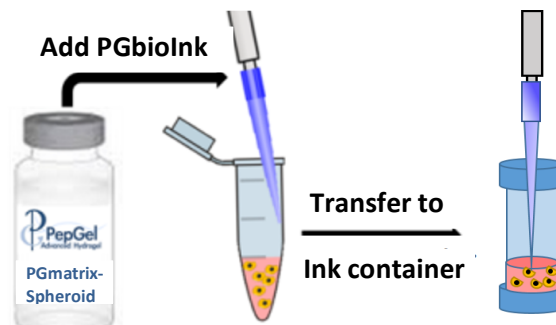

REFERENCE

1. Li Q, Qi G, Liu X, Bai J, Zhao J, Tang G, Zhang YS, Chen-Tsai R, Zhang M, Wang D, Zhang Y. Universal Peptide Hydrogel for Scalable Physiological Formation and Bioprinting of 3D Spheroids from Human Induced Pluripotent Stem Cells. *Advanced Functional Materials*. 2021:2104046.
2. Huang H, Shi J, Laskin J, Liu Z, McVey DS, Sun XS. Design of a shear-thinning recoverable peptide hydrogel from native sequences and application for influenza H1N1 vaccine adjuvant. *Soft Matter*. 2011;7(19):8905-12.
3. Huang H, Ding Y, Sun XS, Nguyen TA. Peptide hydrogelation and cell encapsulation for 3D culture of MCF-7 breast cancer cells. *PloS one*. 2013 Mar 20;8(3):e59482.
4. Li X, Galliher-Beckley A, Huang H, Sun X, Shi J. Peptide nanofiber hydrogel adjuvanted live virus vaccine enhances cross-protective immunity to porcine reproductive and respiratory syndrome virus. *Vaccine*. 2013 Sep 23;31(41):4508-15.
5. Kumar D, Kandl C, Hamilton CD, Shnayder Y, Tsue TT, Kakarala K, Ledgerwood L, Sun XS, Huang HJ, Girod D, Thomas SM. Mitigation of tumor-associated fibroblast-facilitated head and neck cancer progression with anti-hepatocyte growth factor antibody ficlatuzumab. *JAMA Otolaryngology–Head & Neck Surgery*. 2015 Dec 1;141(12):1133-9.
6. Kumar D, Kandl C, Hamilton CD, Shnayder Y, Tsue TT, Kakarala K, Ledgerwood L, Sun X, Huang H, Girod D, Thomas SM. Anti-HGF antibody ficlatuzumab mitigates tumor-associated fibroblast-facilitated head and neck cancer progression. *JAMA Otolaryngology*. 2015;141(12):1133-9.
7. Miller PG, Shuler ML. Design and demonstration of a pumpless 14 compartment microphysiological system. *Biotechnology and bioengineering*. 2016 Oct;113(10):2213-27.
8. Liang J, Sun XS, Yang Z, Cao S. Anticancer drug camptothecin test in 3D hydrogel networks with HeLa cells. *Scientific reports*. 2017 Feb 1;7(1):1-9.
9. Liang J, Liu G, Wang J, Sun XS. Controlled release of BSA-linked cisplatin through a PepGel self-assembling peptide nanofiber hydrogel scaffold. *Amino acids*. 2017 Dec 1;49(12):2015-21.
10. Li L, Liu G, Timashev P, Sun XS, Criswell T, Atala A, Zhang Y. Biofabrication of tissue-specific extracellular matrix proteins to enhance the expansion and differentiation of skeletal muscle progenitor cells. *Applied Physics Reviews*. 2019 Jun 26;6(2):021309.
11. Xu J, Qi G, Sui C, Wang W, Sun X. 3D h9e peptide hydrogel: An advanced three-dimensional cell culture system for anticancer prescreening of chemopreventive phenolic agents. *Toxicology in Vitro*. 2019 Dec 1;61:104599.
12. Thippabhotla S, Zhong C, He M. 3D cell culture stimulates the secretion of in vivo like extracellular vesicles. *Scientific reports*. 2019 Sep 10;9(1):1-4.

13. Zhu Q, Hamilton M, Vasquez B, He M. 3D-printing enabled micro-assembly of a microfluidic electroporation system for 3D tissue engineering. *Lab on a Chip*. 2019;19(14):2362-72.
14. Harrison R, Lockett J, Marsh S, Lugo Leija HA, Salih S, Alkharji R, Sottile V. Magnetically Assisted Control of Stem Cells Applied in 2D, 3D and In Situ Models of Cell Migration. *Molecules*. 2019 Jan;24(8):1563.
15. Carter T, Qi G, Wang W, Nguyen A, Cheng N, Ju YM, Lee SJ, Yoo J, Atala A, Sun XS. Self-Assembling Peptide Solution Accelerates Hemostasis. *Advances in Wound Care*. 2020. (<https://www.liebertpub.com/doi/abs/10.1089/wound.2019.1109>)